I | LABORATORY GLASS BOTTLES AND ACCESSORIES

Due to their high chemical and thermal-shock resistance, DURAN® laboratory bottles are indispensable for routine work in the laboratory. As neutral glass with high hydrolytic resistance, DURAN® belongs to glass type 1 in accordance with EP, JP and USP.

The extensive range of original accessories includes caps for the widest possible range of applications. DURAN® laboratory bottles are completed by suitable pouring rings from different plastics, which enable drip-free working.

Alongside clear glass laboratory bottles, amber (absorbs light up to approx. 500 nm) and plastic-coated (absorbs light up to approx. 380 nm) bottles are also available to protect light-sensitive substances. The PU plastic coating also provides protection against scratching and, in the event of breakage, retains contents and glass splinters.

Usage tips:

- DURAN® laboratory bottles are, with the exception of the pressure-resistant DURAN® pressure plus bottles, not designed for use under pressure or vacuum conditions.
- When sterilising / autoclaving, the screw cap should

only be loosely applied since with a closed bottle pressure equalisation does not take place. An ideal complement is the membrane cap (see pages 20, 34).

- When cleaning in a dishwasher, load such that the glass bottles, especially the threads, are not in contact with each other.
- The printed scale on many items of DURAN® laboratory glassware is indicative with an accuracy of $\pm 10\%$. Therefore the items are not suitable for use as volumetric glassware.
- DURAN® protect bottles can be autoclaved. The maximum permissible temperature is $+135\text{ °C}$ (max. 30 Min.).
- When freezing media, store the bottles in an inclined position and do not fill more than three quarters full. Bear in mind that plastic caps and pouring rings have restricted service temperature ranges.

DURAN® laboratory glass and reagent bottles are provided with a retrace code. Using the eight-character code and the corresponding article number, a batch and quality certificate for the DURAN® laboratory glass bottles can be obtained at www.duran-group.com.

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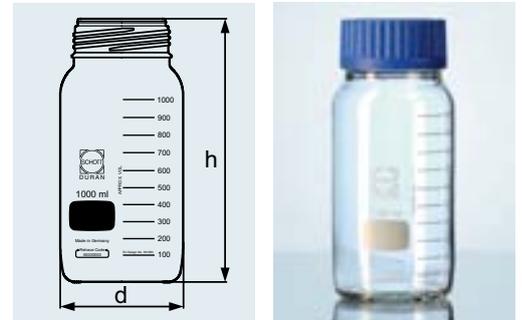
With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue quick release closure (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of closure and pouring ring: +140 °C. Special thread means opening takes less than a turn. The 80 mm wide outer diameter of the bottle mouth permits easy filling and pouring out of powders and viscous substances.

Typical applications: storage, transport, safekeeping and sampling of substances, easy to use with granulated material, powders and viscous media, sampling of hot media.

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DURAN® GLS 80® LABORATORY BOTTLE, WIDE MOUTH
with GLS 80® thread

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DURAN® GLS 80® LABORATORY
BOTTLE, WIDE MOUTH, AMBER
with GLS 80® thread

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue quick release closure (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of closure and pouring ring: +140 °C. Alongside easy handling, UV protection up to 500 nm. Unchanged DURAN® properties within the bottle, as colouration is only to the outer surface. Very uniform, durable and chemically resistant amber colour due to use of innovative technology.

Typical applications: storage, transport and safekeeping of light-sensitive substances, easy to use with granulated material, powders and viscous media.

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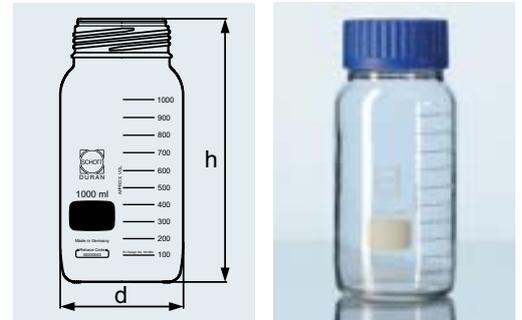
With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue quick release closure (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of closure and pouring ring: +140 °C. Service temperature limit of the PU¹ plastic coating: -30 °C to +135 °C. The coating provides scratch, leak and splinter protection and is ideally suited to both the transport and storage of hazardous media or valuable samples. UV protection up to approx. 380 nm wavelength. High transparency. Suitable for microwaving.

Typical applications: storage, transport and safe handling of hazardous substances. Storage of high value viscous liquids, pastes and powder.

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DURAN® GLS 80® PROTECT
LABORATORY BOTTLE, WIDE MOUTH
with GLS 80® thread, plastic coated

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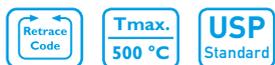
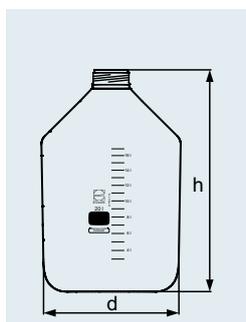
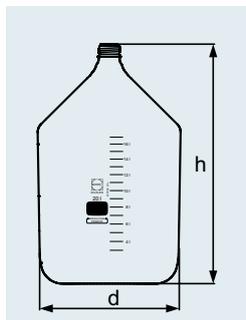
DURAN® PRODUCTION AND STORAGE BOTTLES (CARBOYS)

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These larger sized bottles / carboys are ideal for bulk storage and handling of both liquid and solid intermediates and final formulations. Manufactured from Type I borosilicate 3.3 glass for durable performance and resistance to thermal stress. The glass conforms to American (USP), European (EP) and Japanese (JP) pharmacopoeia standards making the carboys ideal for pharmaceutical production applications. Manufactured with thickened, uniform side walls for higher mechanical strength. Available in two mouth sizes: The versatile DIN standard GL 45 or GLS80® wide mouth for easier cleaning, filling and emptying. Retrace Code for batch traceability and conformance certification. Manufactured from inorganic materials (Certified BST / TSE free). Suitable for high temperature sterilization, depyrogenisation or autoclaving. Feature large, permanent, easy-to-read, white enamel graduations marks. Also available with an external Polyurethane coating for enhanced scratch resistance, and to contain leakage in the event of damage. Available with customized logos, identification labeling or graduations. Supplied without screw cap or pouring ring, but can be used in conjunction caps and connector systems see pages 20 ff and 33 ff for options.

Typical applications: Flat robust base is ideal for mixing processes with large magnetic stir bars.

Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread GL/GLS	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
11 602 00	10 000	45	227,50	410	1
11 601 00	20 000	45	288,50	505	1
11 602 20	10 000	80	227,50	385	1
11 601 10	20 000	80	288,50	480	1

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Permits opening and closing of the DURAN® GLS 80® bottle with only a three-quarter turn. A matching PP¹ pouring ring is also available, permitting clean, drip-free use.

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
QUICK RELEASE CLOSURE FOR DURAN® GLS 80® LABORATORY BOTTLE from PP¹, blue, with lip seal

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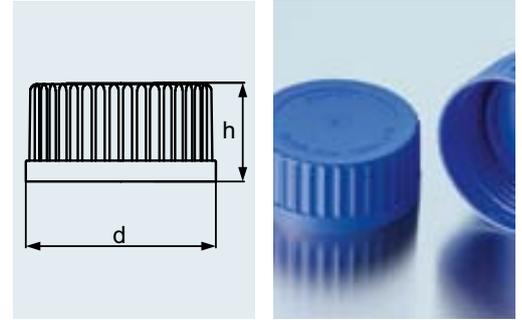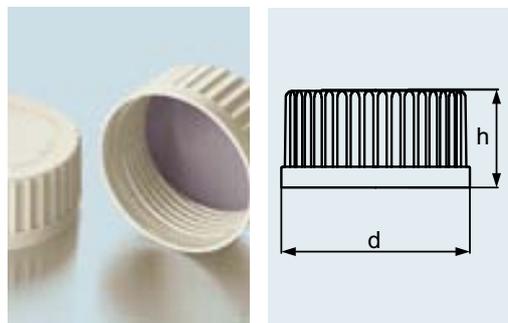
QUICK RELEASE CLOSURE FOR DURAN® GLS 80® LABORATORY BOTTLE
with sealing disc, (PSU compound¹)

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The material used is a special compound based on a polyarylsulphone I. Consequently the chemical, thermal and mechanical properties of the material are noticeably improved and matched to laboratory requirements. Thanks to the thread, the DURAN® GLS 80® bottle can be opened or closed with only a three-quarter turn. The seal, which is coated on both sides with PTFE¹, ensures the bottle can be tightly sealed. A matching PTFE¹ pouring² ring is also available, permitting clean, drip-free use.

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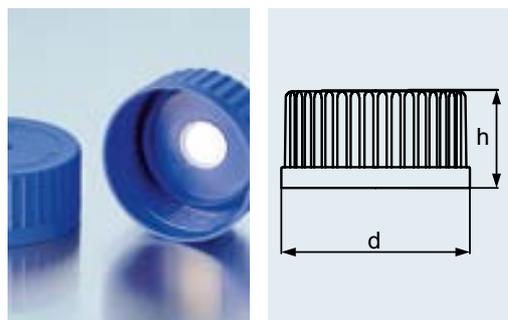
MEMBRANE SCREW CAP FOR DURAN® GLS 80® LABORATORY BOTTLE
from PP¹, blue, with welded-in PTFE¹ membrane
for pressure equalisation

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For GLS 80® thread. Ideal for autoclaving processes because the membrane permits pressure equalisation and screw tight sealing. Hence the risk of contamination is greatly reduced. Ingress of liquids or solids is prevented and the bottle contents remain sterile.

Typical applications: storage or transport of gas generating media, autoclaving of media.

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Materials used: PP¹ and PTFE¹. Flexible modular system. Five different tubing diameters (3.2 mm; 6.0 mm; 8.0 mm; 10.0 mm and 12.0 mm) can be connected. Sterile pressure equalisation is possible by using the syringe filter. Unused ports can be closed with a blind cap.


Typical applications: safe transfer of liquid media within a closed and sterile system (evaporation is reduced).

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GLS 80[®] CONNECTION SYSTEM
screw cap GLS 80[®] with four ports
(GL 18 thread)

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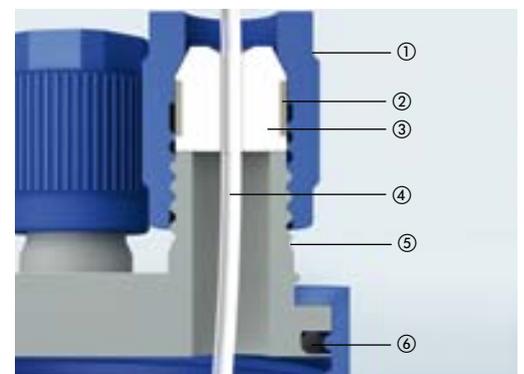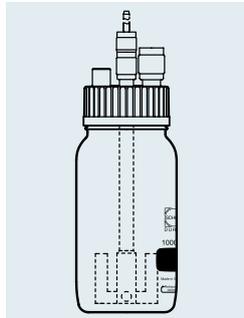
GLS 80® STIRRED REACTOR

materials used: PP'/PTFE'/PEEK'/stainless steel

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The GLS 80® stirred reactor is suitable for a wide range of laboratory mixing processes. The connections (2 x GL 14, 2 x GL 18) provided permit addition or removal of media from the bottle during the mixing process. The whole unit can be autoclaved and is therefore suitable for biological applications. By using components from the GLS 80® connection system, an additional media bottle (OD hose: 1.6 - 12.0 mm)² can be connected or a sterile pressure equalizer attached. Drive for the stirrer is provided by a standard commercial magnetic stirrer. The variable stirrer shaft can be used in DURAN® GLS 80® laboratory glass bottles (1 000 and 2 000 ml) and provides notably improved mixing in comparison with standard magnetic stirring bars. The agitating element is exchangeable and use up to 500 rpm is possible.

Typical applications: mixing of liquids, dissolving of solids, simple fermentation processes.

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DURAN® products with individual labelling

Innovative laser marking, e.g.: Numbering, bar codes, logos or names

No effect on the product performance. More information, see page 229

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² Components for a secure hose connection, see page 21 and 36.

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DURAN® LABORATORY BOTTLE
with DIN thread

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. With proven DURAN® properties. Complete with blue screw cap (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature level of screw cap and pouring ring: +140 °C. For screw cap assortment, see pages 33–42.

Typical applications: storage, sample preparation, transport. Autoclaving media.

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² With specially shaped glass edge for improved pouring out (so that an additional plastic pouring ring is not required).

³ Bottles of 100 – 2 000 ml capacity have a shoulder reinforcement ring corresponding to the bottles nominal capacity.

⁴ Acceptance within ISO 4796 standard has been requested.

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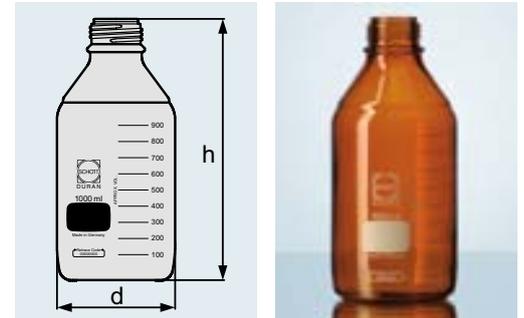
With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. UV protection up to approx. 500 nm wavelength. Unchanged DURAN® properties within the bottle, as colouration is only on the outer surface. Very uniform, durable and chemically resistant amber colour due to use of innovative technology. For screw cap assortment, see pages 33–42.

Typical applications: storage and transport of light-sensitive substances.

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DURAN® LABORATORY BOTTLE, AMBER with DIN thread

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¹ With specially shaped glass edge for improved pouring out (so that an additional plastic pouring ring is not required).

² Bottles of 100 – 2 000 ml capacity have a shoulder reinforcement ring corresponding to the bottles nominal capacity.

³ Acceptance within ISO 4796 standard has been requested.

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DURAN® PROTECT
LABORATORY BOTTLE
 with DIN thread, plastic coated

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Service temperature limit of the PU¹ plastic coating: -30 °C to +135 °C (instructions for autoclaving, see page 217). The coating provides scratch, leak and splinter protection and is ideally suited to both the transport and storage of hazardous media or valuable samples. UV protection up to approx. 380 nm wavelength. High transparency. Suitable for microwaving. Caps and pouring rings see pages 33–42.

Typical applications: storage, transport and safe handling of hazardous or valuable substances.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Ohne Verschluss und Ausgießring, klar Without cap and pouring ring, clear					
10 926 76	25 ²	25	36	70	10
10 926 77	50	32	46	87	10
21 805 24	100 ³	45	56	100	10
21 805 29	150 ⁴	45	62	105	10
21 805 36	250 ³	45	70	138	10
21 805 44	500 ³	45	86	176	10
21 805 51	750 ⁴	45	95	203	10
21 805 54	1 000 ³	45	101	225	10
21 805 63	2 000 ³	45	136	260	10
21 805 69	3 500 ⁴	45	160	295	1
21 805 73	5 000	45	182	330	1

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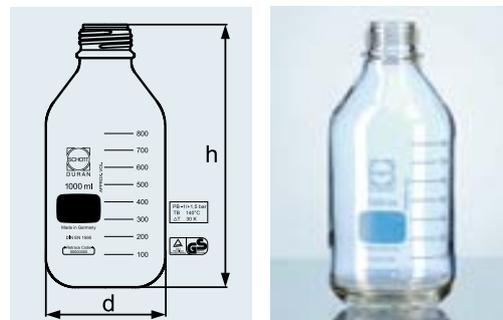
With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable ceramic. Pressure resistance conforms to DIN ISO 1595, confirmed by GS marking (TÜV ID: 0000020716). Vacuum and/or pressure resistant from -1 to +1.5 bar due to modified geometry (based on ISO 4796-1). When pressure loaded the following apply: thermal shock resistance 30 K and maximum usage temperature +140 °C. Blue scale for visual differentiation from the standard laboratory bottle. Also available in amber. See pages 33-42 for spare caps and pouring rings.

Typical applications: safe working under pressure or vacuum, sampling under pressure, storage of gas generating media.

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DURAN® PRESSURE PLUS LABORATORY BOTTLE¹
with DIN thread, GL 45

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Ohne Verschluss und Ausgießring, klar Without cap and pouring ring, clear					
10 922 34	250	45	70	138	10
10 922 35	500	45	86	176	10
21 810 54	1 000	45	101	225	10
Braun Amber					
10 943 67	250	45	70	138	10
10 943 68	500	45	86	176	10
21 816 54	1 000	45	101	225	10

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¹ Bottle with plastic coating is available on demand.

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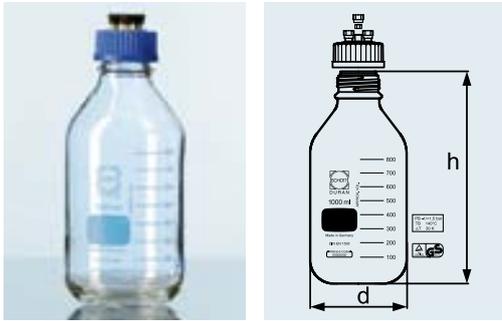
DURAN® HPLC BOTTLE
with DIN thread, GL 45

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable ceramic. Complete system comprising: DURAN® pressure plus laboratory bottle with 4-port screw cap (PP¹); four screw connections (black, M8 thread) and silicone seals. Connection of different hose diameters (1.6 mm and 3.2 mm) as well as sterile pressure equalisation sets (syringe filter 0.2 µm) is possible. Unused ports can be sealed with silicone blanking seals.

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³ Replacement bottle, see page 27.

⁴ Including silicone seals.

⁵ GL 45 screw cap, 4 x M8 screw cap, 12 x silicone seal.

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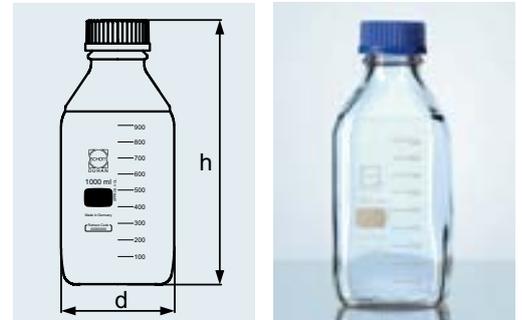
With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue screw cap (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of screw cap and pouring ring: +140 °C. Ergonomic handling due to angular shape, highly stable, good stackability. Alongside proven DURAN® properties, a space saving of 44% in comparison with standard laboratory bottles (example applies to 100 ml bottles). Screw caps are also available in the following colours: green, yellow and grey (see pages 33–42).

Typical applications: space-saving storage, space-saving transport.

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**DURAN® LABORATORY BOTTLE,
SQUARE
with DIN thread**

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with pouring ring and cap from TpCh260¹ (similar to PFA). The premium cap with its PTFE¹ coated silicone seal is colourless and temperature resistant from –196 °C to +260 °C. Together with proven DURAN® properties, TÜV tested thermal shock resistance of 160 K, confirmed by GS-marking (TÜV ID: 0000020715). USP/FDA conformity for the entire system comprised of bottle, screw cap and pouring ring. DMF available upon request (DMF no. 19757). Accurate scale: ± 5%. Additional graduations as well as additional opposing scale simplify reading off.

Typical applications: Due to its properties, ideal for applications in the pharmaceutical industry, handling of aggressive media, sterilisation processes (hot air and dry sterilisation) and depyrogenation.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
11 270 75	100	45	56	105	10
11 270 76	250	45	70	143	10
11 270 77	500	45	86	181	10
11 270 78	1 000	45	101	230	10

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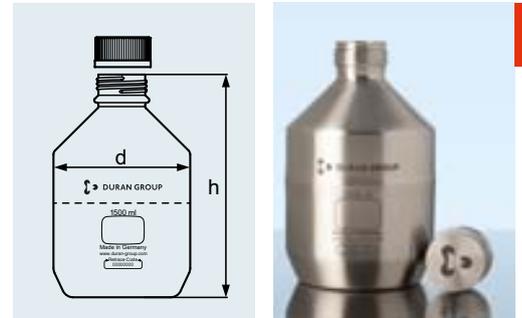
The unbreakable DURAN Group steel bottle is ideal for storage applications where glass is not applicable, due to the risk of breakage or chemical incompatibility. Manufactured from corrosion resistant AISI Type 316L (1.4404) stainless steel, and hygienically constructed with all welded seams. The bottle has a smooth inner surface finish that corresponds to IIIc (DIN 17441) with rounded inner edges for ease of cleaning. Polished and brushed durable exterior finish. The bottle has a GL 45 thread with a built-in pouring lip. Supplied without a cap, stainless steel cap is available. It is fully compatible with all the GL 45 caps (see page 33 ff).

Typical applications: Hazardous materials laboratory container for storage of liquids, intermediates, or solid product. Storage of precious materials, such as high purity fine chemicals, pharmaceutical or cosmetic products. Storage of light sensitive materials.

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DURAN GROUP STAINLESS STEEL LABORATORY BOTTLE
with DIN thread, GL 45


Best.-Nr. Cat. No.		Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungs- einheit/ Stück Pack/Quantity
29 901 60 0	DURAN Group GL 45 Edelstahlflaschen DURAN Group GL 45 Stainless Steel Laboratory Bottle	1 500	45	120	201	1
29 911 28	DG Edelstahlverschluss (316L), mit platinver- netzter Silikondichtung und PTFE Beschichtung, GL 45 Stainless steel (316L) cap, with PTFE surfaced Platinum-catalyzed silicon liner, GL 45		45	50	27	1

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DURAN GROUP STAINLESS STEEL SHIPPING BOTTLE
with DIN thread GL 45, UN certified

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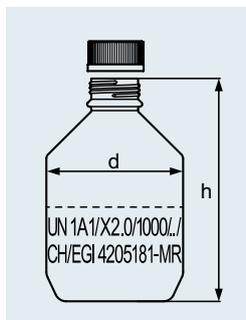The unbreakable DURAN Group steel bottle is ideal for the storage and safe shipping of dangerous liquid goods such as solvents, and reagents, without the need for additional protective outer packaging¹. Manufactured from corrosion resistant AISI Type 316L (1.4404) stainless steel, and hygienically constructed with all welded seams. The bottle has a smooth inner surface finish that corresponds to IIIc (DIN 17441) with rounded inner edges for ease of cleaning. Polished and brushed durable exterior finish. The bottle has a GL 45 thread with a built-in pouring lip. Supplied complete with GL 45 Stainless Steel screw cap and PTFE surfaced Platinum-catalyzed silicon liner, and UN certification number.

Typical applications: Container for the shipping of hazardous or non-hazardous liquids. Storage or shipping of precious liquids, such as high purity fine chemicals, pharmaceutical or cosmetic products. Storage of light sensitive liquids.

Best.-Nr. Cat. No.		Inhalt Capacity ml	Verpackungs- einheit/ Stück Pack/Quantity
29 901 60 5	DURAN Group GL 45 (Transport-) Edelstahlflasche, UN zertifiziert ¹ , komplett DURAN Stainless Steel GL 45 Shipping Bottle, UN certified ¹ , complete	1 500	1

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¹ Certified to UN standards for the carriage of liquids classified as dangerous goods in UN packing Groups II (medium danger) and III (low danger). It is suitable for liquids with a relative density of 2.0 or less. International regulations are subject to change, it is the user's responsibility for complying with all applicable laws and regulations.

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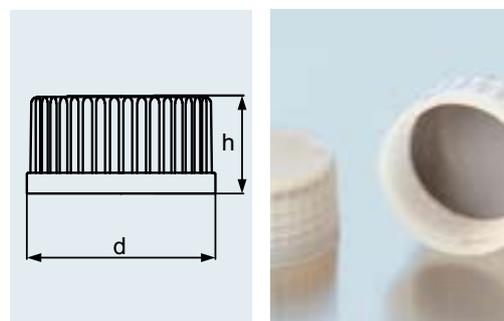
Cap is moulded from a dye-free polymer to avoid risk of leaching. Wide usable temperature range from -196 °C to $+260\text{ °C}$. Very high chemical resistance. Complete with colourless PTFE¹ coated silicone sealing disk for high leak tightness. Registered as Type III Drug Master File in conjunction with Premium bottle (DMF no. 19757). Copy available on request. The matching TpCh260¹ pouring ring permits drip-free pouring is available separately. Replacement PTFE¹ coated silicone sealing disks² are available.

Typical applications: Due to its exceptional properties, it is ideal for applications in the pharmaceutical industry, storage of aggressive materials, and demanding sterilisation processes such as hot air sterilisation or depyrogenation.

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PREMIUM CAP
from TpCh260¹ (similar to PFA) with
PTFE¹ coated silicone seal²

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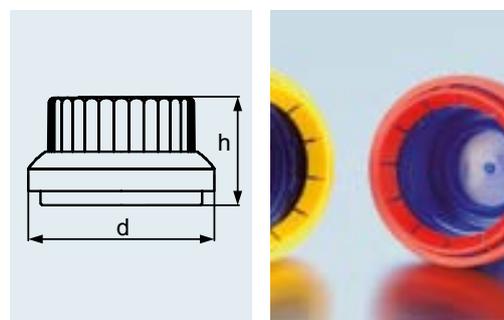
The tamper-evident screw cap is available with lip seal or PTFE¹ coated silicone seal². The integral plastic ring tears when the cap is first opened and is retained on the bottle neck. Therefore it can be clearly seen whether the bottle retains its original seal. After initial removal, can be used as “normal” screw cap.

Typical applications: secure storage and transport/shipping of valuable media.

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TAMPER-EVIDENT SCREW CAP
from PP¹, for laboratory bottles with DIN thread

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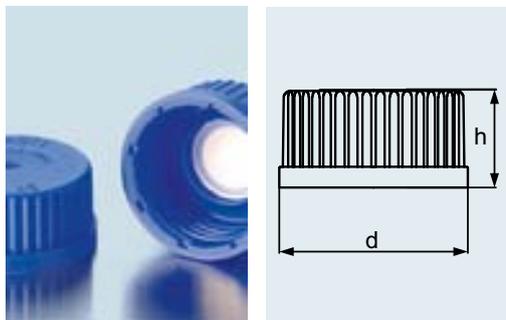
MEMBRANE SCREW CAP
 from PP¹, blue with welded-in
 PTFE¹ membrane for pressure equalisation

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*For all GL 45, GL 32 and GL 25 threads.
 Ideal for autoclaving processes because
 the membrane permits pressure equalisation
 and tight sealing, greatly reducing the risk of
 contamination. Ingress of liquids or solids is
 prevented and the bottle contents remain
 sterile.*

*Typical applications: storage or transport
 of gas generating media, autoclaving of
 media.*

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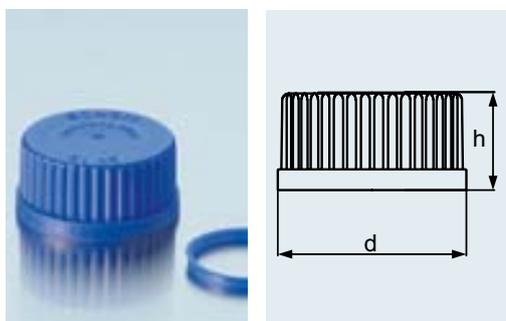
*Available in the colours blue, green, yellow
 and grey with matching colour pouring
 rings. Distinguishing media types is simplified
 and interchanging of screw caps and carry
 over of substances is practically eliminated.*

*Typical applications: visual aid to the safe
 identification of different media.*

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
High leak tightness through use of PTFE¹ coated silicone seal³. More chemically resistant than PP¹ screw cap. A matching ETFE¹ pouring ring is also available, permitting clean, drop-free use.

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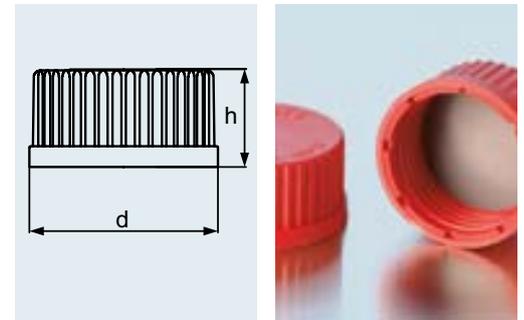¹ For chemical and thermal resistance, see page 222.

² For screw connection caps with aperture, silicone seals with aperture and septa, see pages 179–181.

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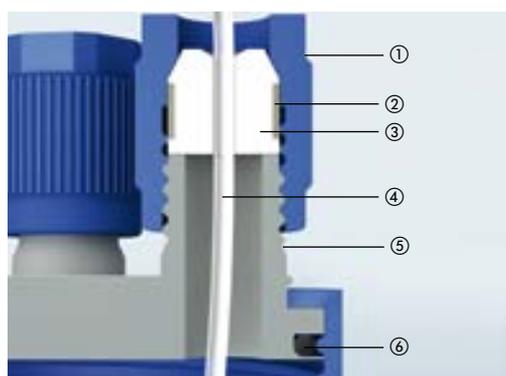
CONNECTION SYSTEM

screw cap GL 45 with two or three ports (GL 14 thread)

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Materials used: PP¹ and PTFE¹. Flexible modular system. Four different tubing diameters (1.6 mm; 3.0 mm; 3.2 mm and 6.0 mm) can be connected. Sterile pressure equalisation is possible through use of syringe filter. Unused ports can be provided with a blind cap.

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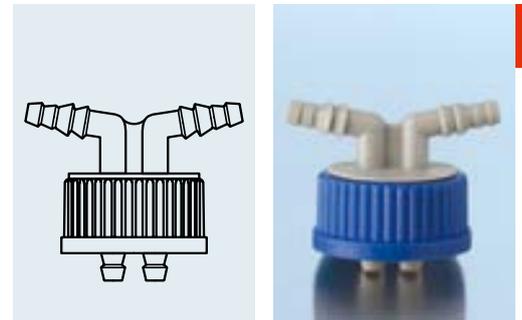
GL45 cap connection system for the easy connection of flexible tubing to the DURAN® GL45 bottles. The two polypropylene connectors have angled top and straight underside connectors. The cap is ideal for use with soft elastic tubing that has an inner diameter of 6 – 9 mm (such as Viton®, Tygon®, or other silicone tubing). An optional venting connector is available with, or without a syringe filter.

Highly versatile as the screw cap is based on the standard GL 45 thread. Grey polypropylene cap center rotates freely, allowing the bottle to be unscrewed without the need to disconnect the tubing. The angled top connectors prevent kinks in the tubing. Temperature resistant up to +140 °C. Fully autoclavable and washer-proof.


Typical applications: Possible biotech applications include the transfer of sterile media from one container to another using a peristaltic pump.

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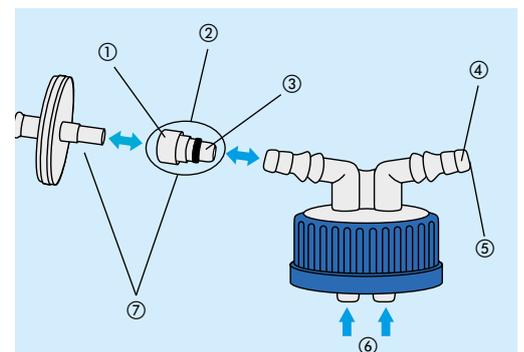verwendete Materialien PP¹/PTFE¹/PEEK¹/
 Edelstahl

GL 45 STIRRED REACTOR

materials used PP¹/PTFE¹/PEEK¹/stainless steel

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Self contained stirring system ideal for mixing processes. Stirrer shaft can be adjusted to the optimal position in either a 500 or 1,000 ml DURAN® GL 45 bottles. Drive with standard magnetic stirres up to 500 rpm. Significantly improved mixing performance compared to standard magnetic stirring bars. Multi-connector cap is compatible with the proven DURAN® connection systems; tubing with 1.6 mm to 6.0 mm bores can be used to add or remove liquid or gas. Fully autoclavable. Parts in contact with media conform to FDA requirements. Available separately or with bottle.

Typical applications: mixing chemicals, media or cultures in the DURAN® GL 45 laboratory bottle.

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DURAN Group Bottle-top dispensers make the serial dispensing of liquids directly from DURAN® laboratory glass bottles safer and quicker.


Reliable and reproducible liquid dispensing due to a very precise borosilicate 3.3 glass cylinder. Volume settings can be quickly and accurately adjusted.

Simple dispensing operation, even with one hand. Dispenser rotates freely 360° on the bottle to a convenient position. Materials of construction offer good chemical resistance. Adaptable to different bottles threads (adaptors for GL 32, GL 38, S 40 or GL 28, GL 45, S 40) and bottle heights with a telescopic intake tube. PFA sealing lip prevents crystallization build-up. The modular design is easy to dismantle and clean.

Supplied with certificate of performance, and multi-lingual instruction manual.

Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread GL	Adapter Adaptor	Verpackungseinheit/ Stück Pack/Quantity
29 115 03	0,25 – 2,5	32	GL 28, GL 45, S40	1
29 115 06	0,5 – 5,0	32	GL 28, GL 45, S40	1
29 115 09	1,0 – 10,0	32	GL 28, GL 45, S40	1
29 115 14	2,5 – 25,0	45	GL 32, GL 38, S40	1
29 115 17	5,0 – 50,0	45	GL 32, GL 38, S40	1
29 115 24	10,0 – 100,0	45	GL 32, GL 38, S40	1

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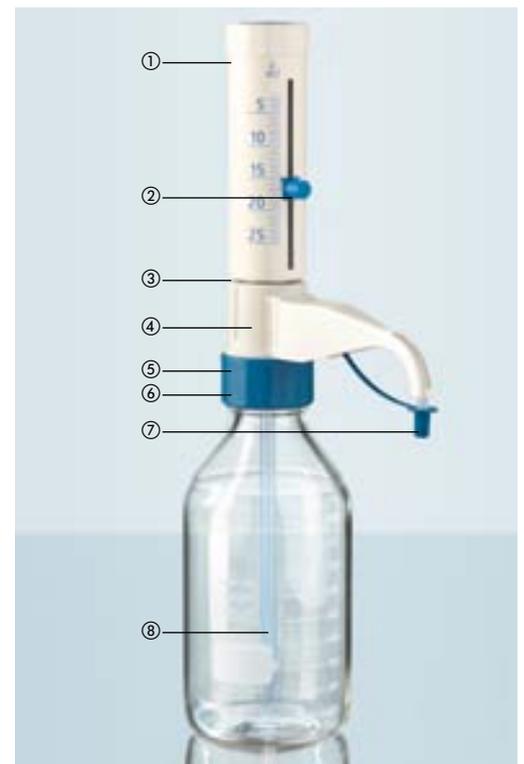DG SAFETY CAP
screw cap GL 45 with one, two or three ports

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High-quality and flexible connection system, specifically designed for handling aggressive, gas-generating media. Materials used: PP¹, PBT¹ and PTFE¹. All parts coming into contact with media are from PTFE¹. Three different hose diameters (1.6 mm; 2.3 mm; 3.2 mm) can be connected. Pressure equalisation is possible through use of membrane filter. Outgassing of media is reduced by use of a one-way valve that decreases media loss and protects the user from vapours. A free moving PTFE¹ core prevents rotation of the hoses when screwing onto the DURAN[®] bottle. Unused ports can be provided with a blind cap. Optionally available with stopcock to prevent hoses running dry or air intakes during bottle changeover. DG safety caps are also available for bottles with standard ground joint.

Typical applications: safe transfer of aggressive media within a closed system (evaporation is reduced).

Best.-Nr. Cat. No.	Bezeichnung Description	DIN-Gewinde DIN Thread GL	Verpackungseinheit/ Stück Pack/Quantity
11 684 01 ²	DG Safety Cap I (1 Anschluss) DG Safety Cap I (1 port)	45	1
11 684 02 ²	DG Safety Cap II (2 Anschlüsse) DG Safety Cap II (2 ports)	45	1
11 684 13 ²	DG Safety Cap III (3 Anschlüsse) DG Safety Cap III (3 ports)	45	1
11 684 14 ²	DG Safety Cap I (1 Anschluss), mit Absperrhahn DG Safety Cap I (1 port), with stopcock	45	1
11 684 15 ²	DG Safety Cap II (2 Anschlüsse), mit Absperrhähnen DG Safety Cap II (2 ports), with stopcocks	45	1
11 684 16 ²	DG Safety Cap III (3 Anschlüsse), mit Absperrhähnen DG Safety Cap III (3 ports), with stopcocks	45	1
11 684 17 ³	DG Safety Cap für NS 29/32 DG Safety Cap for NS 29/32		1

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¹ For chemical and thermal resistance, see page 222.

² Complete with GL 45 screw cap, connection screws, conical nipple (ID 3.2 mm, PTFE), ventilation valve and membrane filter.

³ Complete with screw cap for NS 29/32, connection screws, conical nipple (ID 3.2 mm, PTFE), ventilation valve and membrane filter.

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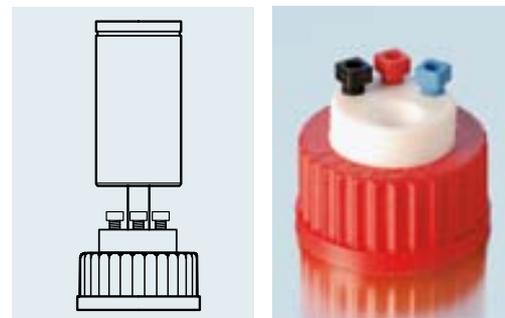
The activated carbon filter prevents the escaping of hazardous gases and thus increases safety in the laboratory. Materials used: PBT¹ and PTFE¹. Optional special connection for installations with a leakage system. A free moving PTFE¹ core prevents rotation of the hoses when screwing onto the DURAN® bottle. Unused ports can be provided with a blind cap.

Typical applications: safe storage and collection of liquid laboratory waste.

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DG SAFETY WASTE CAP
additional connection option for activated carbon filter and leak hose, GL 45

Best.-Nr. Cat. No.	Bezeichnung Description	DIN-Gewinde DIN Thread GL	Verpackungseinheit/ Stück Pack/Quantity
11 684 18 ²	DG Safety Waste Cap DG Safety Waste Cap	45	1
11 684 20 ³	DG Safety Waste Cap für Schlauchverbindung DG Safety Waste Cap for hose connector	45	1

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¹ For chemical and thermal resistance, see page 222.

² Complete with GL 45 screw cap, connection screws and conical nipple (ID 3.2 mm, PTFE).

³ Complete with GL 45 screw cap, connection screws, conical nipple (ID 3.2 mm, PTFE) and connection for leak hose.

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
*Matching cap, see pages 44–45.
A maximum usage temperature of +100 °C is recommended. Thermal shock resistance 30 K. Hydrolytic class 3.*

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SCREW CAP BOTTLE, SQUARE, FROM SODA-LIME GLASS
narrow neck with thread

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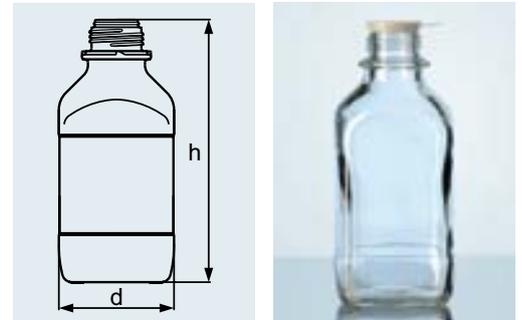
*Without pouring ring and dust cover cap¹.
A maximum usage temperature of +100 °C is recommended. Thermal shock resistance 30 K. Hydrolytic class 3.*

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SCREW CAP BOTTLE, SQUARE, FROM SODA-LIME GLASS
wide neck with thread

Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Klartglas <i>Clear glass</i>					
23 820 17	50	32 ²	48	70	10
23 820 24	100	32 ²	49	111	10
23 820 36	250	45 ²	64	146	10
23 820 44	500	54 ²	76	173	10
23 820 54	1 000	60 ²	97	213	10
Braunglas <i>Amber glass</i>					
23 826 24	100	32 ²	49	111	10
23 826 36	250	45 ²	64	146	10
23 826 44	500	54 ²	76	173	10
23 826 54	1 000	54 ²	97	213	10

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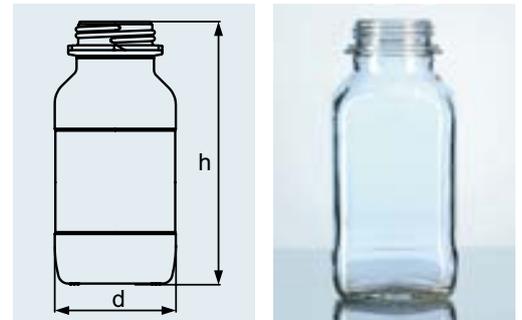SCREW CAP BOTTLE, ROUND, AMBER
FROM SODA-LIME GLASS
 with thread

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Complete with pouring ring and dust cover
 cap. A maximum usage temperature of
 +100 °C is recommended. Thermal shock
 resistance 30 K. Hydrolytic class 3.

Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Braunglas Amber glass					
23 835 66 5	2 500	45 ¹	139	283	1

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TAMPER-EVIDENT SCREW CAP
 from PP¹, blue, for soda-lime square screw
 cap bottle, narrow neck

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SECURITY SCREW CAP
 from PP¹ with valve², red, for soda-lime
 square screw cap bottles, narrow neck

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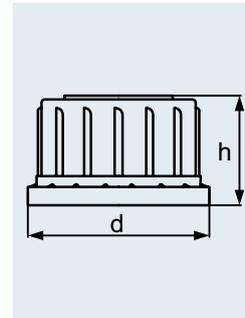
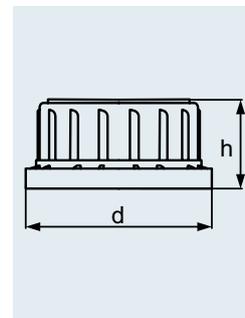
TAMPER-EVIDENT SCREW CAP
 from PP¹, blue, for soda-lime
 square screw cap bottles, wide neck

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With precision ground neck. All glass components, therefore also suitable for storage of aggressive media, which could attack plastic parts.

Typical application: storage of powders.

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DURAN® REAGENT BOTTLE, WIDE NECK

neck with standard ground joint,
with standard ground glass flat-head stopper

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With precision ground neck. All glass components, therefore also suitable for storage of aggressive media, which could attack plastic parts. Also available in amber. Unchanged DURAN® properties within the bottle, as colouration is only to the outer surface. Very uniform amber colour thanks to innovative technology, moreover long durability and high chemical resistance of the amber colouring.

Typical applications: storage of powders.

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¹ Narrow neck bottles with standard ground joint, 1 000 ml, also suitable for use as burette bottles, see page 124.

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With precision ground neck. All glass components, therefore also suitable for storage of aggressive media, which could attack plastic parts. Also available in amber. Unchanged DURAN® properties within the bottle, as colouration is only on the outer surface. Very uniform, durable and chemically resistant amber colour due to use of innovative technology.

Typical application: storage of liquids.

Best.-Nr. Cat. No. klar/clear	Best.-Nr. Cat. No. braun/amber	Inhalt Capacity ml	d mm	h mm	Hals Neck NS	Verpackungseinheit/ Stück Pack/Quantity
Hals mit Normschliff und NS-Glasstopfen Neck with standard ground joint and glass flat-head stopper						
21 165 08 ¹		10	28	52	10/19	10
21 165 14	21 168 14	25	36	64	12/21	10
21 165 17	21 168 17	50	42	80	14/15	10
21 165 24	21 168 24	100	52	96	14/15	10
21 165 36	21 168 36	250	70	130	19/26	10
21 165 44	21 168 44	500	86	164	24/29	10
21 165 54	21 168 54	1 000	107	200	29/32	10
21 165 63	21 168 63	2 000	134	248	29/32	10
21 165 73	21 168 73	5 000	182	323	45/40	1
21 165 86	21 168 86	10 000	227	398	60/46	1
21 165 91	21 168 91	20 000	288	492	60/46	1

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

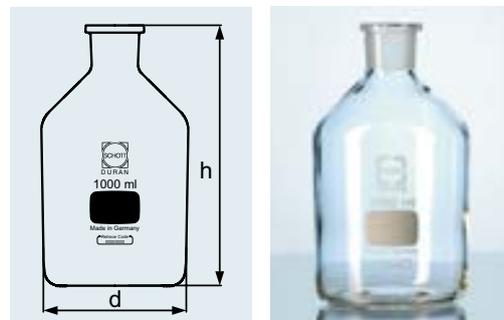
DURAN® REAGENT BOTTLE, NARROW NECK

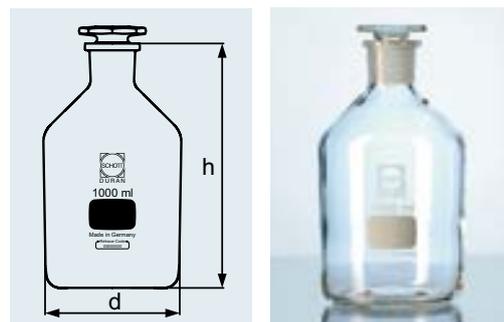neck with standard ground joint, with standard ground glass flat-head stopper

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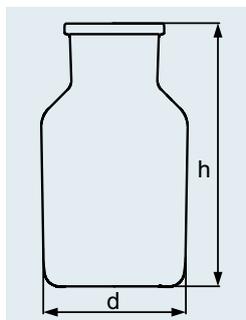REAGENT BOTTLE, WIDE NECK
 FROM SODA-LIME GLASS
 neck with standard ground joint

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Without stopper¹. A maximum usage temperature of +100 °C is recommended. Thermal shock resistance 30 K. Hydrolytic class 3.

Best.-Nr. Cat. No. klar/clear	Best.-Nr. Cat. No. braun/amber	Inhalt Capacity ml	d mm	h mm	Hals Neck NS	Verpackungseinheit/ Stück Pack/Quantity
Hals mit Normschliff, ohne Stopfen Neck with standard ground joint, without stoppers						
23 184 17		50	44	79	24/20	10
23 184 24	23 187 24	100	52	97	29/22	10
23 184 36	23 187 36	250	71	129	34/24	10
23 184 44	23 187 44 ²	500	86	164	45/40	10
23 184 54	23 187 54	1 000	107	200	60/46	10

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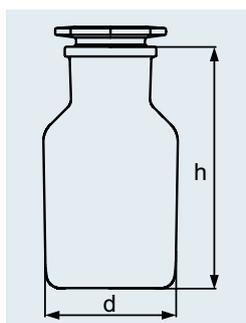
Note on using DURAN® glass stoppers with soda-lime glass bottles: If the bottle and the stopper have a temperature difference greater than 30 °C, the glass stoppers can become stuck!

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REAGENT BOTTLE, WIDE NECK
 FROM SODA-LIME GLASS
 neck with standard ground joint,
 with standard ground glass flat-head stopper¹

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A maximum usage temperature of +100 °C is recommended. Thermal shock resistance 30 K. Hydrolytic class 3.

Best.-Nr. Cat. No. klar/clear	Best.-Nr. Cat. No. braun/amber	Inhalt Capacity ml	d mm	h mm	Hals Neck NS	Verpackungseinheit/ Stück Pack/Quantity
Hals mit Normschliff und NS-Glasstopfen Neck with standard ground joint and glass flat-head stopper						
23 185 17		50	44	79	24/20	10
23 185 24	23 188 24	100	52	97	29/22	10
23 185 36	23 188 36	250	71	129	34/24	10
23 185 44	23 188 44 ²	500	86	164	45/40	10
23 185 54	23 188 54	1 000	107	200	60/46	10

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REAGENT BOTTLE, NARROW NECK
FROM SODA-LIME GLASS
neck with standard ground joint


Best.-Nr. Cat. No. klar/clear	Best.-Nr. Cat. No. braun/amber	Inhalt Capacity ml	d mm	h mm	Hals Neck NS	Verpackungseinheit/ Stück Pack/Quantity
Hals mit Normschliff Neck with standard ground joint						
	23 167 17	50	42	80	14/15	10
23 164 24	23 167 24	100	52	96	14/15	10
23 164 36	23 167 36	250	72	130	19/26	10
23 164 44	23 167 44	500	89	165	24/29	10
23 164 54 ¹	23 167 54 ¹	1 000	110	200	29/32	10

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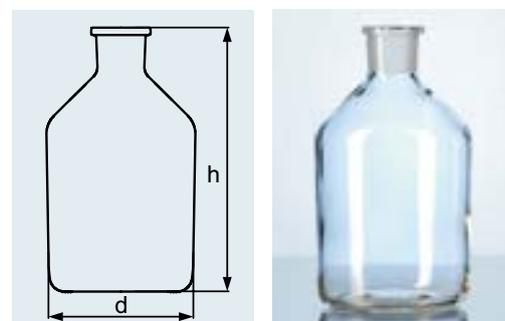
REAGENT BOTTLE, NARROW NECK
FROM SODA-LIME GLASS
neck with standard ground joint,
with standard ground glass flat-head stopper

Best.-Nr. Cat. No. klar/clear	Best.-Nr. Cat. No. braun/amber	Inhalt Capacity ml	d mm	h mm	Hals Neck NS	Verpackungseinheit/ Stück Pack/Quantity
Hals mit Normschliff und NS-Glasstopfen Neck with standard ground joint and glass flat-head stopper						
	23 168 17	50	42	80	14/15	10
23 165 24	23 168 24	100	52	96	14/15	10
23 165 36	23 168 36	250	72	130	19/26	10
23 165 44	23 168 44	500	89	165	24/29	10
23 165 54	23 168 54	1 000	110	200	29/32	10

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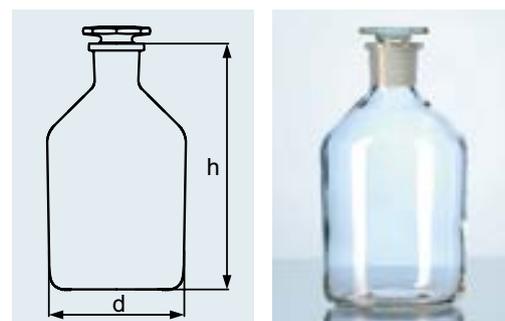GLASS STOPPER, AMBER
with standard ground joint, octagonal

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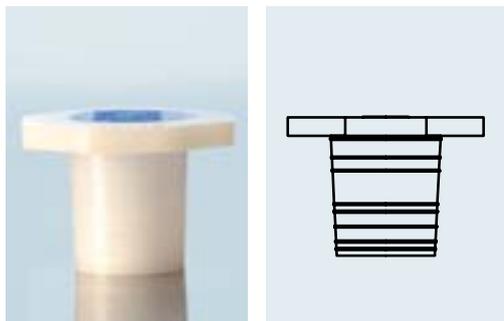
PLASTIC STOPPER
from polyethylene¹

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GLASS STOPPER
with standard ground joint, octagonal

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GLASS STOPPER
with short ground joint, octagonal

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DURAN® GLASS STOPPER
ground conical, for reagent bottles
(oxygen bottles according to Winkler)

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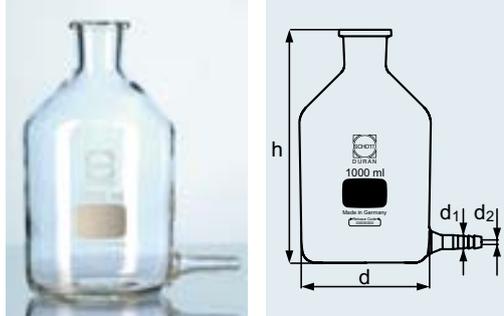
Note on using DURAN® glass stoppers with **soda-lime** glass bottles: If the bottle and the stopper have a temperature difference greater than 30 °C, the glass stoppers can become stuck!

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DURAN® ASPIRATOR (LEVELLING)
BOTTLE
with plain neck and bottom sidearm

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
Useful as a delivery or storage container for solutions. Tubular sidearm (Diameter: Outer d_1 : 11 mm, Inner d_2 : 5 mm). Outlet facilitates attachment of flexible tubing.

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
24 708 36	250	73	130	10
24 708 44	500	89	164	10
24 708 54	1 000	111	200	1

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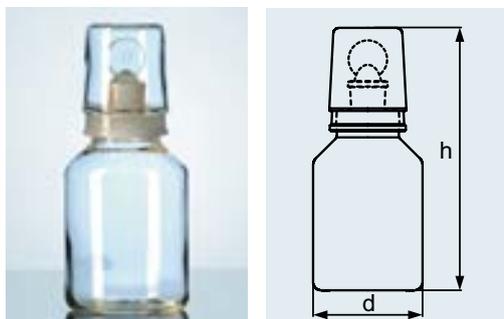DURAN® ACID BOTTLE¹
with standard ground "pennyhead" stopper,
conical shoulders, interchangeable glass cap

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In addition to the ground stopper, a glass cap with ground joint is supplied. This provides an improved seal and protection against acid vapours.

Best.-Nr. Cat. No.	Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	NS	Verpackungseinheit/ Stück Pack/Quantity
klar clear	braun amber					
21 275 24	21 275 24 6	100	55	145	19/17	10
21 275 36	21 275 36 6	250	75	180	19/17	10
21 275 44	21 275 44 6	500	82	220	24/20	10
21 275 54	21 275 54 6	1 000	109	260	29/32	10

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DROPPING BOTTLE
FROM SODA-LIME GLASS
with interchangeable clear glass pipette¹ standard
ground joint, complete with rubber teats²

Best.-Nr. Cat. No.	Best.-Nr. Cat. No.	Inhalt Capacity ml	h mm	NS	Verpackungseinheit/ Stück Pack/Quantity
klar clear	braun amber				
23 270 17	23 270 17 6	50	79	14/15	10
23 270 24	23 270 24 6 ³	100	105	14/15	10

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
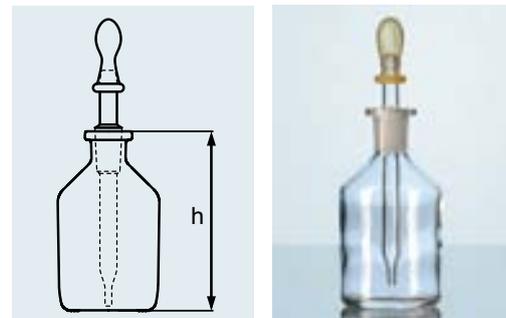
Dosing of liquids is possible via an outlet.

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DURAN® ASPIRATOR (LEVELLING)
BOTTLE
tubulated with standard ground joint,
without stoppers, neck unground


Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Bodentubus Tubulature NS	Verpackungseinheit/ Stück Pack/Quantity
24 701 44	500	86	164	19/26	10
24 701 54	1 000	107	200	19/26	10
24 701 63	2 000	134	249	19/26	10
24 701 73	5 000	182	320	29/32	1
24 701 86	10 000	228	398	29/32	1
24 701 91	20 000	289	492	29/32	1

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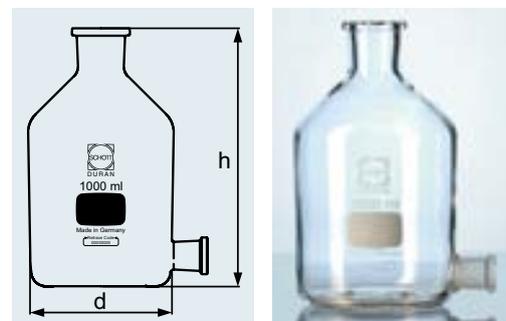DURAN® ASPIRATOR (LEVELLING)
BOTTLE

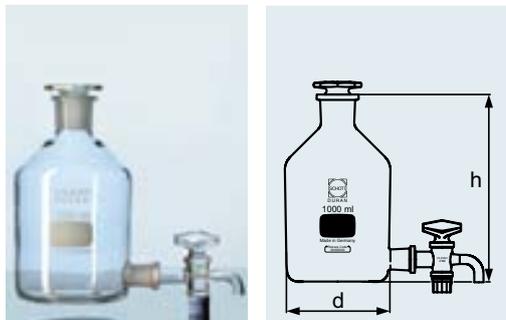*tubulated with standard ground joint,
 complete with standard ground stopcock
 and standard ground stopper*

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Dosing of liquids is possible via a stopcock.

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Bodentubus Tubulature NS	Hals Neck NS	Orig. Pack./Stück Stand. pack/quantity
24 702 44	500	86	164	19/26	24/29	10
24 702 54	1 000	107	200	19/26	29/32	10
24 702 63	2 000	134	249	19/26	29/32	10
24 702 73	5 000	182	320	29/32	45/40	1
24 702 86	10 000	228	398	29/32	60/46	1
24 702 91	20 000	289	492	29/32	60/46	1

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STOPCOCK WITH STANDARD
GROUND JOINT FOR ASPIRATOR
BOTTLE

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DURAN® ASPIRATOR (LEVELLING)
BOTTLE

*neck with DIN thread GL 45,
 tubulator with GL 32*

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*Complete with screw connection cap¹,
 silicone seal and stopcock with PTFE²
 spindle. Dosing of liquids is possible via a
 stopcock.*

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
24 703 54	1 000	45	101	225	1
24 703 63	2 000	45	136	260	1
24 703 73	5 000	45	182	330	1
24 703 86	10 000	45	230	410	1

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² For chemical and thermal resistance, see page 222.


Best.-Nr. Cat. No.	Bohrung Hole mm	für Inhalt for Capacity ml	Orig. Pack./Stück Stand. pack/quantity
24 147 03	6	1 000 + 2 000	1
24 147 04	8	5 000 + 10 000	1

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STOPCOCK FOR ASPIRATOR BOTTLE
with PTFE¹ spindle, for GL 32 screw thread

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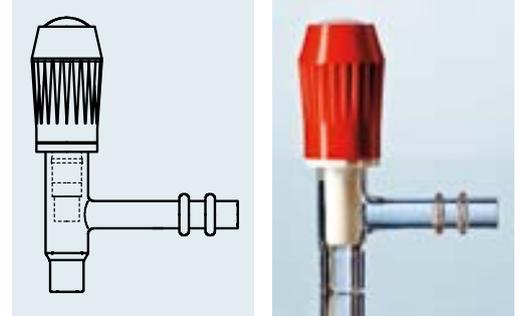PTFE-SILICONE¹ SEAL
for GL 32 screw thread

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